



Grindometers – Fineness of Grind Gauges VF2110, VF2111, VF2112, VF2113

Manual

1 PRODUCT DESCRIPTION

Precision instrument to determine particle size and fineness of many materials like paints, lacquers, pigments, filler, chocolate etc.. Most TQC Grindometers have double grooves with graded slopes graduated in three different parameters: Microns, NS (Hegman) and PCU (North). Gauge and beveled scraper are made of hardened stainless steel and have an accuracy of 2 µm.

2 SPECIFICATIONS AND ARTICLE NUMBERS

The base: 175 x 60 x 12mm with 120mm groove length

Tolerance: ±2µm Stainless steel

The scraper: 75 x 38 x 8mm shaped to be hold easily.

Stainless steel





VF2110 Grindometer FM15/2 Din-Iso Range: 0-15 micron, 10-8,5 PCU, 8-6,8 Hegman Graduation: 1,5 micron Double groove

VF2111 Grindometer FM25/2 Din-Iso Range: 0-25 Micron, 10-7,5 PCU, (north), 8-6 Hegman (NS) Graduation: 2,5 micron Double groove

VF2112 Grindometer FM50/2 Din-Iso Range: 0-50 Micron, 10-5 PCU, (north), 8-4 Hegman (NS) Graduation: 5 micron Double groove VF2113 Grindometer FM100/2 Din-Iso

Range: 0-100 Micron, 10-0 PCU, (north), 8-0 Hegman (NS)

Graduation: 10 micron Double groove

VF2108 Grindometer FM100/1 Din-lso Range: 0-100 Micron, 8-0 Hegman (NS) Graduation: 10 micron

Single groove

3 STANDARDS

ASTM D 1210, ASTM D 1316, JIS K 5600-2-5, ISO 1524, DIN EN 21524, BS 3900-C6

teprüfung fineriess of grind madflijhneid Mantfelnheit drying fine droogtijd Trockenzeit abrasion slijtvastheid Abrieblestigkeit inspection accessories t lanz & 'Haze' corrosion carrosie Karrosion coating thickness gauges coatingdikte meters Schichtdicken messgeräte flash point vlampunt Flammpun the testapparatuur Klimaprüfung appearance uiterlijk Erscheinungsbild inspection kits inspectie kits Inspektionskoffer elasticity & resistance deformat ances weegschalen Waagen surface cleanliness oppervlaktezuiverheid Oberflächenreinheit film application filmapplicatie Filmaufziehgeräte coat

 TQC bv.
 2761 JP Zevenhuizen
 Tel. +31 (0)180-631344
 E-mail
 info@tqc.eu

 Nijverheidscentrum 14
 The Netherlands
 Fax +31 (0)180-632917
 http
 www.tqc.eu

vf2110-vf2113_m44.doc - Page 2 of 2

4 TAKING A READING

When using the gauge, take care of not to damage the surface of the gauge or the edges of the scraper.

Ensure the surface of gauge and edge of the scraper is clean from material residue, oil, etc.

Perform a preliminary test to determine the size of gauge most suitable for the fineness of grind characteristics of the material being tested.

- 1. Place the gauge on a flat, horizontal and non-slip surface, with the zero mark on the scale closest to the user.
- 2. Place a suitable amount of the material in the deep end of each groove.
- 3. Place the scraper on the surface of the gauge behind the material. Use both hands to hold the scraper as shown in **Picture 1.**
- 4. Pull the scraper along the length of the gauge at a constant speed and apply sufficient downward Pressure to clean excess material from the edges of the gauge. This operation takes approx. 1 to 2 seconds.
- 5. View the drawn out material within the next 3 seconds. This avoids inaccurate testing due to evaporation of the material. The material should be viewed at right angles to the length of the groove and at an angle of 20° to 30° with the surface of the gauge.
- 6. Find a band across the grooves of 3mm wide which contains 5 to 10 particles of the material as shown in **Picture 2.** Read the position of the upper limit of this band on the scale and record this value.
- 7. Use a suitable solvent to clean the gauge and also the scraper.
- 8. Perform 2 more tests and calculate the average value of the results. The average value is the fineness of grind of the material.



Picture 1.



Picture 2.

5 MAINTENANCE

- Always clean the instrument after use with a suitable solvent.
- Never clean the instrument by any mechanical means such as a wire brush or abrasive paper. This may
 cause, just like the use of aggressive cleaning agents, permanent damage.
- The instruments have to be protected from rust when it is not in use. Rust can appear on the instrument when it is used only occasionally and when it is been handled by a user with sweaty hands.
- Always dry the instrument and scraper after use to protect against rust, and apply a thin layer of oil to the surface of the instrument and scraper before storage.
- Always store the instrument in its pouch when not in use.
- Check regularly whether the gauge and the scraper are worn or damaged.

6DISCLAIMER

The information given in this manual is not intended to be exhaustive and any person using the product for any purpose other than that specifically recommended in this manual without first obtaining written confirmation from us as to the suitability of the product for the intended purpose does so at his own risk. Whilst we endeavour to ensure that all advice we give about the product (whether in this manual or otherwise) is correct we have no control over either the quality or condition of the product or the many factors affecting the use and application of the product. Therefore, unless we specifically agree in writing to do so, we do not accept any liability whatsoever or howsoever arising for the performance of the product or for any loss or damage (other than death or personal injury resulting from our negligence) arising out of the use of the product. The information contained in this manual is liable to modification from time to time in the light of experience and our policy of continuous product development.

teprüfung ihreress of grind waalfijnneid Mantfelnheit dryng Time aroogtlijd Trackerzelt abrasion slijfvasineid Abriebtestigkeit inspection accessories lanz & 'Haze' corrosion corrosie Korrosion coating thickness gauges coatingdikte meters Schichtdicken messgeräte flash point vlampunt Flammpur the testapparatuur Klimaprüfung appearance uiterlijk Erscheinungsbild inspection kits inspectie kits Inspektionskoffer elasticity & resistance deforma ances weegschalen Waagen surface cleanliness oppervlaktezuiverheid Oberflächenreinheit film application filmapplicatie Filmaufziehgeräte coa