Phosphate CHEMets® Kit


K-8530/R-8515: 2 - 30 ppm **K-8515/R-8515**: 10 - 150 ppm


Safety Information

Read MSDS (available at www.chemetrics.com) before performing this test procedure. Wear safety glasses and protective gloves.


Test Procedure

- 1. Fill the sample cup to the 25 mL mark with the sample to be tested (fig 1).
- Place the CHEMet ampoule, tip first, into the sample cup. Snap the tip. The ampoule will fill leaving a bubble for mixing (fig 2).
- To mix the ampoule, invert it several times, allowing the bubble to travel from end to end
- Dry the ampoule and wait 5 minutes for color development.
- 5. Obtain a test result using the comparator.
 - a. For K-8530 (fig 3): Place the ampoule, flat end first, into the comparator. Hold the comparator up toward a source of light and view from the bottom. Rotate the comparator until the best color match is found.
 - b. For K-8515 (fig 4): Place the ampoule between the color standards until the best color match is found.

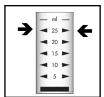

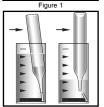



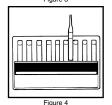

Test Method

The Phosphate CHEMets®1 test method employs the vanadomolybdophosphoric acid chemistry.^{2,3} In an acidic solution, ortho-phosphate reacts with ammonium molybdate and ammonium vanadate to produce a yellow colored complex in direct proportion to the phosphate concentration.

Condensed phosphates (pyro-, meta-, and other polyphosphates) and organically bound phosphates do not respond to this test. Sulfide, thiosulfate, and thiocyanate will cause low test results.

- CHEMets is a registered trademark of CHEMetrics, Inc. U.S. Patent No. 3.634.038
- 2. APHA Standard Methods, 21st ed., method 4500-P C (2005)
- 3. ASTM D 515 82, Phosphorus in Water, Test Method C

Visit www.chemetrics.com to view product demonstration videos.

Always follow the test procedure above to perform a test.



www.chemetrics.com 4295 Catlett Road, Midland, VA 22728 U.S.A. Phone: (800) 356-3072; Fax: (540) 788-4856 E-Mail: orders@chemetrics.com

Jan. 13, Rev. 5