Compact vertical table-top-type

Compact vertical table-top-type with high ability

Designed for the purpose of "One set for one person." Compact body of W37 × D30 × H30 cm. And yet, capacity of three 1,000 ml vial bottles or four 500 ml vial bottles (In case of using wire basket, two or three). High ability of reliable sterilization of max. 127°C.

Up-opening lid & front operation design occupies minimum space

The lid opens and closes lightly with one hand, yet securely due to the self-closing silicon rubber packing. The up-opening lid and only front operation design, realize ideal space saving. You can reasonably lay out your laboratory.

Operation-start is one-touch of the start button

By just pushing the start button, you get fully automatic operation from start to finish. The same parts as middle or large models are used the automatic air exhaust device guarantees perfect sterilization under pure steam exposure, and the exact pressure/temperature regulator. The timer begins to move only after the set pressure/temperature is reached, and cuts off electricity after the set time.

One touch setting of sterilizing temperature

Before operation, you can select and set sterilizing temperature by the control knob. So you need not adjust it in sterilizing operation. Also change of temperature (ex. 121 ← 110°C) is one-touch.

Prevention of water lack

The sensitive thermostat mounted on the surface of immersion-type sheathed heater, cuts off immediately the electrical source in case of water lack, and the buzzer sounds continuously.



Safety

Shiny polished SUS-304 stainless steel chamber and double lid with pressed thick plate arm, passed through water pressure test (3.4 kg/cm².... twice as high pressure as safety valve).

Pressure safety valve with enough inner diameter and stainless spring, operates surely at 1.7 kg/cm2 in case of abnormal pressure rise.

Safety breaker (main switch) cuts off the electrical source in case of short circuit or over current of the heater, etc.


Receiver tank of exhaust drain

This receiver tank is set inside the front body, so there is no trouble of falling down. Drained water in the tank is easily poured out to a sink, because it can be pulled out.





My autoclave usable any time

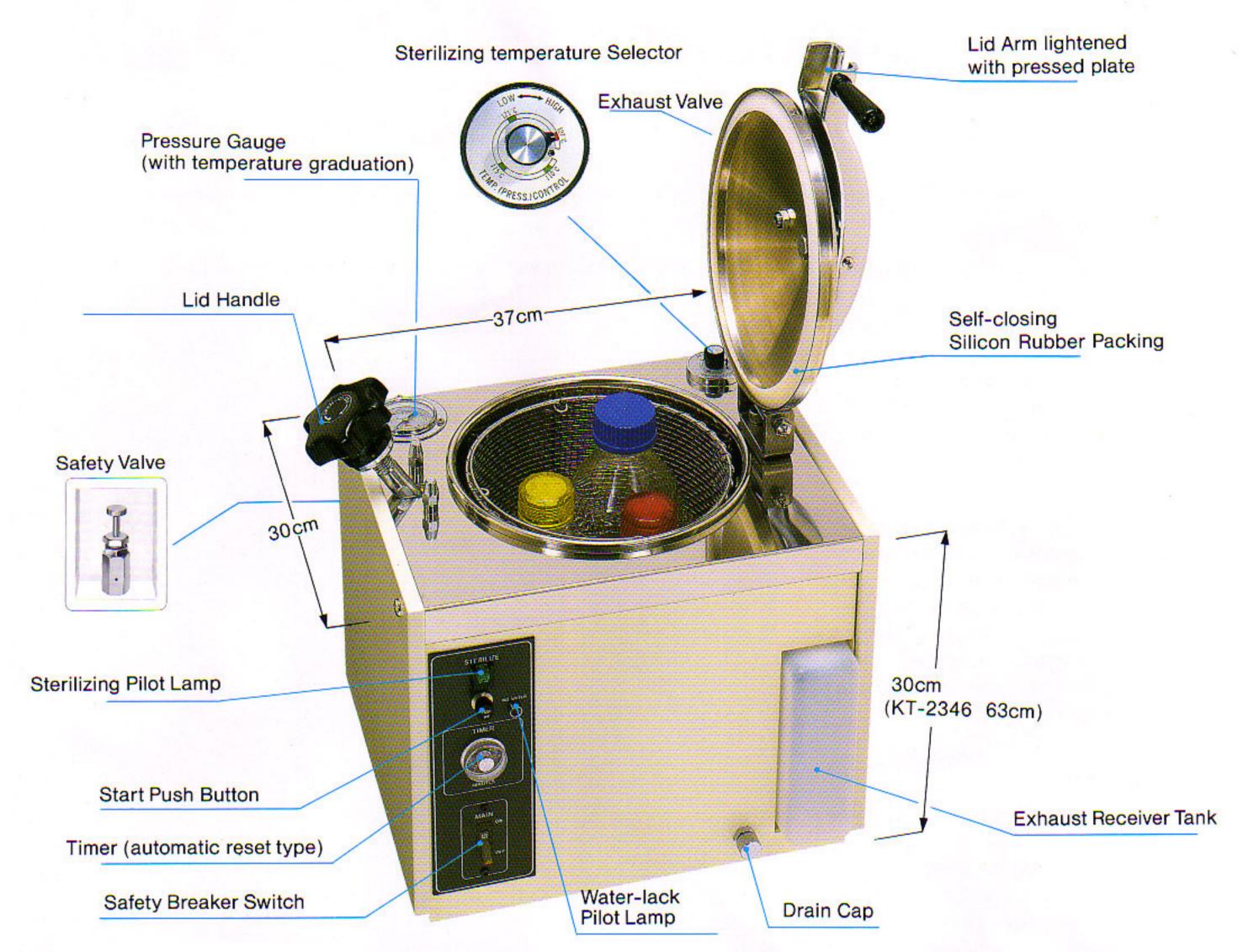

Spec	ific	atic	n
------	------	------	---

	Model	KT-2322	KT-2346
	Chamber	23 cm Ø × 22 cm	23 cm Ø × 46 cm
(Inner dia × usable depth)		(10 <i>l</i>)	(20 <i>l</i>)
	Temperature range	110 ~ 127°C	
ability	One touch set temp.	110, 115, 121, 127°C	
	Usable press. range	0.2 ~ 1.6 kg/cm ² G	
	Max. pressure	1.8 kg/cm ² G	
	Operation start	Push button switch	
system	Timer	60 minutes (Auto. reset type)	
	Press. regulator	Precise press. switch	
	Air exhaust device	Heat operating valve	
	Temp. selector	Dial system	
	Exhaust valve	Screw valve with teflon sheet	
	Drain valve	Cap with silicon sheet	
material	Chamber & lid	SUS304 stainless steel	
	Lid packing	Silicon rubber	
nate	Lid handle	Heat-resisting plastic	
-	Outer plate	Steel with baked enamel finish	

Electricity Outer size (cm) Weight Stainless basket	AC100 ~ 240 \ W37 × D30 × H30 16 kg 21cmø×20cm, 1pc	V, 1∅, 50/60 Hz W37 × D30 × H73 24 kg 21cm∮×40cm, 1 pc.	
Outer size (cm)	$W37 \times D30 \times H30$	W37 × D30 × H73	
		Charles of the Control of the Control	
Electricity	AC100 ~ 240 \	√, 1∅, 50/60 Hz	
THE REPORT OF THE PARTY OF THE	AC100 ~ 240 V, 1Ø, 50/60 Hz		
Heater	1.2 kW		
ain receiver tank	1 <i>l</i>		
standard	by the Ministry of Labor		
Pressure vessel	[Summary pressure vessel] in Japanese standard authorized		
Water test press.	3.4 kg/cm ² G		
Salety valve	(Stainless spring system)		
Safety valve	1.7 kg/cm ² G		
Electric breaker	(Slight leakage breaker: option)		
Water lack preventor	Over current breakter		
	(Heater cut off, Buzzer & Red lamp)		
	by a sensitive thermostat and relay		
	preventor Electric breaker Safety valve Water test press. Pressure vessel standard ain receiver tank	by a sensitive the (Heater cut off, But Over curre (Slight leakage 1.7 kg (Stainless sp. Water test press. Pressure vessel standard Prescription of the distribution of the standard of the cut off, But Over curre (Slight leakage 1.7 kg (Stainless sp. Stainless sp. Stainless sp. Standard of the cut off, But Over curre (Slight leakage 1.7 kg (Stainless sp. Stainless sp. Stainless sp. Standard of the cut off, But Over curre (Slight leakage 1.7 kg (Stainless sp. Stainless sp. Stainless sp. Stainless sp. Standard of the cut off, But Over curre (Slight leakage 1.7 kg (Stainless sp. Stainless sp. Stainl	

Accessories

 Stainless wire basket 		
	KT-2322 21φ×20 cm 1 pc KT-2346 21φ×40 cm 1 pc	
Stainless buck	et	
	(option)	
21 cm Ø × 21	cm	

Dressing drum



(option)

Please specify the voltage when you order.



ALP Co., Ltd.

3-3-10,Midorigaoka Hamura-shi Tokyo,Japan TEL.81-42-579-0531 FAX.81-42-579-0533