

Methods

Iron is present in nature in the form of its oxides, or in combination with silicon or sulfur. The soluble iron content of surface waters rarely exceeds 1 mg/L, while ground waters often contain higher concentrations. The National Secondary Drinking Water Standard for iron is 0.3 mg/L, as iron concentrations in excess of 0.3 mg/L impart a foul taste and cause staining. High concentrations in surface waters can indicate the presence of industrial effluents or runoff.

Iron contamination in oil field brines are typically a result of corrosion processes of iron-containing metallic components and equipment. Accumulation of insoluble iron salts in a brine completion fluid can result in substantial formation damage and can significantly affect the productivity of an oil well. Quantifying total iron in brine is critical.

The Phenanthroline Method (total & soluble; total & ferrous)

References: APHA Standard Methods, 21st ed., Method 3500-Fe B (2005). ASTM D 1068-77, Iron in Water, Test Method A.

With the Phenanthroline Method, ferrous iron reacts with 1,10-phenanthroline to form an orange-colored chelate. To determine total iron, thioglycolic acid solution is added to reduce ferric iron to the ferrous state. The reagent formulation minimizes interferences from various metals. Results are expressed as ppm (mg/L) Fe.

The PDTS Method (total)

Reference: J. A. Tetlow and A. L. Wilson, "Determination of Iron in Boiler Feedwater," *Analyst*, 1958.

CHEMetrics' colorimetric method for determining total iron uses thioglycolic acid to dissolve particulate iron and to reduce iron from the ferric to the ferrous state. Ferrous iron then reacts with PDTS (3-(2-pyridyl)-5,6-bis(4-phenylsulfonic acid)-1,2,4-triazine disodium salt) in acid solution to form a purple-colored chelate. Results are expressed as ppm (mg/L) Fe.

The Ferric Thiocyanate (Iron in Brine)

Reference: D. F. Boltz and J. A. Howell, eds., *Colorimetric Determination of Nonmetals*, 2nd ed., Vol. 8, p. 304 (1978).

The Iron in Brine test employs the ferric thiocyanate chemistry. In an acidic solution, hydrogen peroxide oxidizes ferrous iron. The resulting ferric iron reacts with ammonium thiocyanate forming a red-orange colored

thiocyanate complex, in direct proportion to the iron concentration.

The method expresses total iron content in units of mg/L. By simply dividing the measured mg/L by the density of the brine expressed in units of kg/L, the ppm value can be obtained in mg/kg.

Visual Kits

Range: 0-1 & 1-10 ppm MDL: 0.05 ppm / Method: Phenanthroline	
Iron (total & ferrous) CHEMets Kit	Cat# K-6210
CHEMets Refill, 30 ampoules	R-6201
Activator Solution Pack, six 10 mL bottles	A-6000 ¹
Low Range Comparator 0, 0.1, 0.2, 0.3, 0.4, 0.6, 0.8, 1.0 ppm	C-6001
High Range Comparator 1, 2, 3, 4, 5, 6, 7, 8, 10 ppm	C-6010
Kit comes in a plastic case and contains everything needed to perform 30 tests: Refill, Low and High Range Comparators, Activator Solution, 25 mL sample cup, instructions, and MSDS.	

Range: 0-1 & 1-10 ppm MDL: 0.05 ppm / Method: Phenanthroline	
Iron (total & soluble) CHEMets Kit	Cat# K-6010
CHEMets Refill, 30 ampoules	R-6001
Activator Solution Pack, six 10 mL bottles	A-6000 ¹
Low Range Comparator 0, 0.1, 0.2, 0.3, 0.4, 0.6, 0.8, 1.0 ppm	C-6001
High Range Comparator 1, 2, 3, 4, 5, 6, 7, 8, 10 ppm	C-6010
Kit comes in a plastic case and contains everything needed to perform 30 tests: Refill, Low and High Range Comparators, Activator Solution, 25 mL sample cup, instructions, and MSDS.	

Range: 0-30 & 30-300 ppm MDL: 5 ppm / Method: Phenanthroline	
Iron (total & soluble) VACUettes Kit	Cat# K-6010D
VACUettes Refill, 30 ampoules	R-6001D
Activator Solution Pack, six 10 mL bottles	A-6000 ¹
Low Range Comparator 0, 5, 7.5, 10, 15, 20, 25, 30 ppm	C-6001D
High Range Comparator 30, 60, 90, 120, 150, 175, 200, 250, 300 ppm	C-6010D
Kit comes in a plastic case and contains everything needed to perform 30 tests (except distilled water): Refill, Low and High Range Comparators, Activator Solution, dilutor snapper cup, micro test tube, 5 mL sample cup and top, instructions, and MSDS.	

Range: 0-60 & 60-600 ppm MDL: 10 ppm / Method: Phenanthroline	
Iron (total & soluble) VACUettes Kit	Cat# K-6010A
VACUettes Refill, 30 ampoules	R-6001A
Activator Solution Pack, six 10 mL bottles	A-6000 ¹
Low Range Comparator 0, 10, 15, 20, 30, 40, 50, 60 ppm	C-6001A
High Range Comparator 60, 120, 180, 240, 300, 350, 400, 500, 600 ppm	C-6010A
Kit comes in a plastic case and contains everything needed to perform 30 tests (except distilled water): Refill, Low and High Range Comparators, Activator Solution, dilutor snapper cup, micro test tube, 5 mL sample cup and top, instructions, and MSDS.	

Range: 0-1200 & 1200-12,000 ppm MDL: 200 ppm / Method: Phenanthroline	
Iron (total & soluble) VACUettes Kit	Cat# K-6010C
VACUettes Refill, 30 ampoules	R-6001C
Activator Solution Pack, six 10 mL bottles	A-6000 ¹
Low Range Comparator 0, 200, 300, 400, 600, 800, 1000, 1200 ppm	C-6001C
High Range Comparator 1200, 2400, 3600, 4800, 6000, 7000, 8000, 10,000, 12,000 ppm	C-6010C
Kit comes in a plastic case and contains everything needed to perform 30 tests (except distilled water): Refill, Low and High Range Comparators, Activator Solution, dilutor snapper cup, micro test tube, 5 mL sample cup and top, instructions, and MSDS.	

Range: 0-120 & 120-1200 ppm MDL: 20 ppm / Method: Phenanthroline	
Iron (total & soluble) VACUettes Kit	Cat# K-6010B
VACUettes Refill, 30 ampoules	R-6001B
Activator Solution Pack, six 10 mL bottles	A-6000 ¹
Low Range Comparator 0, 20, 30, 40, 60, 80, 100, 120 ppm	C-6001B
High Range Comparator 120, 240, 360, 480, 600, 700, 800, 1000, 1200 ppm	C-6010B
Kit comes in a plastic case and contains everything needed to perform 30 tests (except distilled water): Refill, Low and High Range Comparators, Activator Solution, dilutor snapper cup, micro test tube, 5 mL sample cup and top, instructions, and MSDS.	

Range: 0-100 & 100-1000 mg/L MDL: 5 mg/L / Method: Ferric Thiocyanate	
Iron in Brine CHEMets Kit	Cat# K-6002
CHEMets Refill, 30 ampoules	R-6002
Acidifier Solution Pack, six 20 mL bottles	A-6001 ¹
Activator Solution Pack, six 20 mL bottles	A-6002 ¹
Low Range Comparator 0, 10, 20, 30, 40, 60, 80, 100 mg/L	C-6002
High Range Comparator 100, 200, 300, 400, 500, 600, 700, 800, 1000 mg/L	C-6012
Kit comes in a plastic case and contains everything needed to perform 30 tests (except distilled water): Refill, Low and High Range Comparators, Acidifier Solution, Activator Solution, 50 mL sample cup with cap, 1.0 mL syringe (2 ea), instructions, and MSDS.	

If no shelf-life is listed for a product, then the shelf-life is at least 2 years.





Instrumental Kits

V-2000 Multi-Analyte Photometer

(See page 12 for instrumental features)

Range: 0.10-2.50 ppm

Method: PDTS

Iron (total) Vacu-vials Kit

Cat#

K-6023

Kit comes in a cardboard box and contains everything needed to perform 30 tests: thirty ampoules, Activator Solution, 25 mL sample cup, ampoule blank, instructions, calibration table, and MSDS.

Range: 0.20-6.00 ppm

Method: Phenanthroline

Iron (total & ferrous) Vacu-vials Kit

Cat#

K-6203

Kit comes in a cardboard box and contains everything needed to perform 30 tests: thirty ampoules, Activator Solution, 25 mL sample cup, ampoule blank, instructions, calibration table, and MSDS.

Range: 0.20-6.00 ppm

Method: Phenanthroline

Iron (total & soluble) Vacu-vials Kit

Cat#

K-6003

Kit comes in a cardboard box and contains everything needed to perform 30 tests: thirty ampoules, Activator Solution, 25 mL sample cup, ampoule blank, instructions, calibration table, and MSDS.

Range: 1.0-25.0 ppm

Method: Phenanthroline

Iron (total & soluble) Vacu-vials Kit

Cat#

K-6013

Kit comes in a cardboard box and contains everything needed to perform 30 tests: thirty ampoules, Activator Solution, 25 mL sample cup, ampoule blank, instructions, calibration table, and MSDS.

Vacu-vials Kits require the use of the V-2000 Photometer or a spectrophotometer capable of accepting a 13 mm diameter round cell. Instrument sold separately.

Kit Components common to Iron

Description	Cat#
Sample Cup Pack, 25 mL (6 ea)	A-0013
Micro Test Tube Pack (10 ea)	A-0015
Dilutor Snapper Cup Pack (6 ea)	A-0018
Ampoule Blank Pack (5 ea)	A-0023
Syringe Pack, 1.0 mL (6 ea)	A-0027
Sample Cup & Cap Pack, 50 mL (6 ea)	A-0058
Sample Cup & Top Pack, 5 mL (6 ea)	A-0105

¹The accessory pack supplies enough solution to perform at least 100 CHEMet or Vacu-vial tests and 42 VACUette tests. A-6000 Activator Solution is required for total iron analysis only.

Instructions are posted on our website.

